



Configuring the NI 6216-BNC USB DAQ Box for use with MiDAS DA

Date Published: January 2013


Abstract

This application note describes the procedures for installing jumper cables and labels on a model 6216-BNC USB DAQ box from National Instruments for use with MiDAS DA.

Instructions

MIDAS DA is designed to work with a wide range of hardware from National Instruments. In order to properly route synchronization and trigger signals to the appropriate BNC connectors on hardware not purchased from Xcitex, users must attach jumpers to the screw terminal block.


1. Re-label the BNC connector on the top right named USER to **Sync Out**, as shown below.

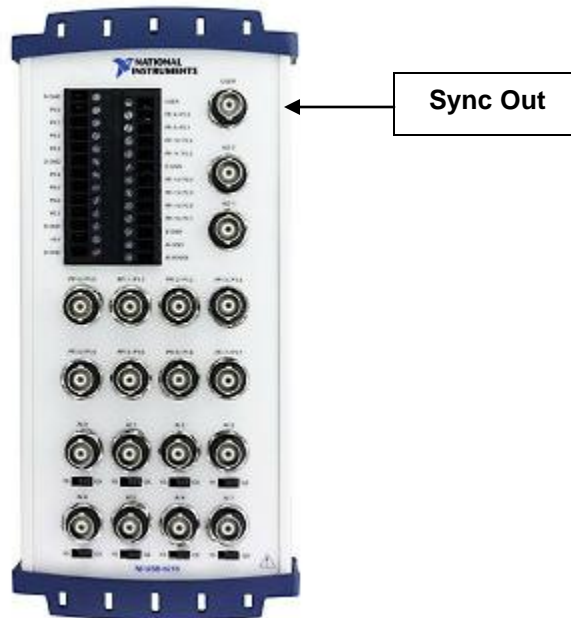


- Using a BNC to 2-wire adapter (BNC pigtail) as shown below,




connect the black lead (-) to **D GND** and connect the red lead (+) to **PFI 11 / P2.3**, as shown in the following pictures. Label this new BNC connector **Trigger In**.

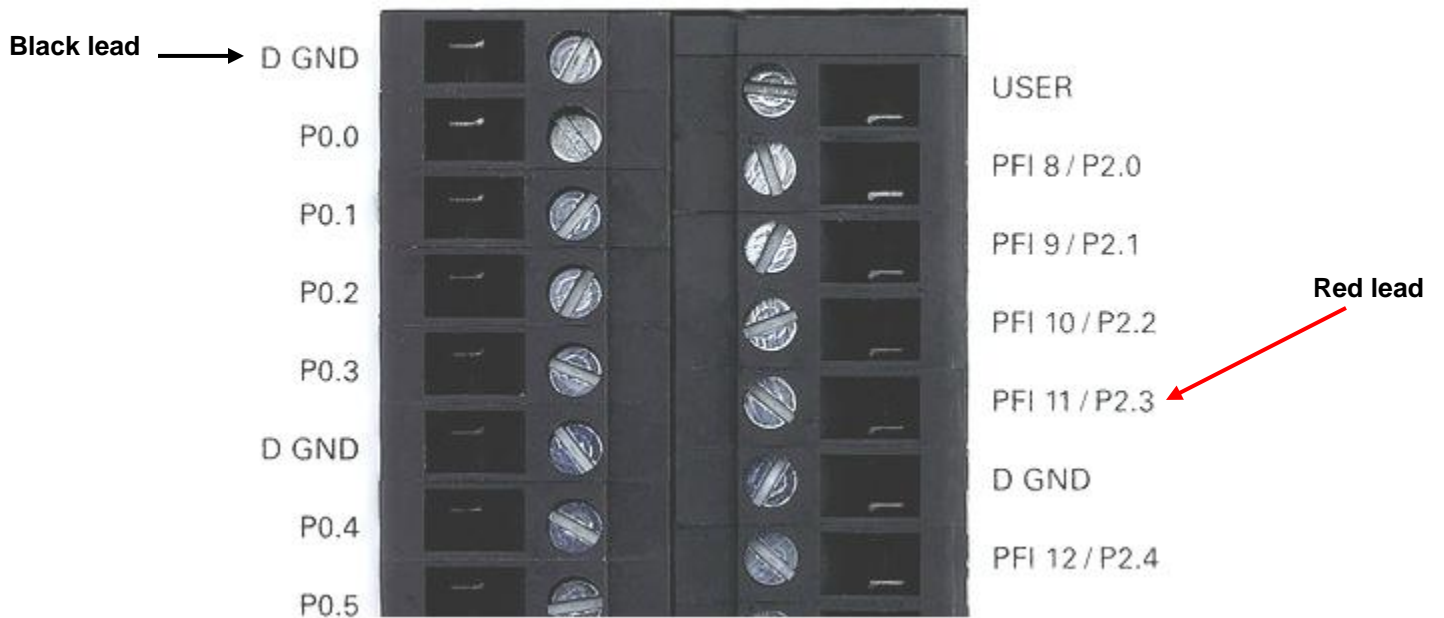


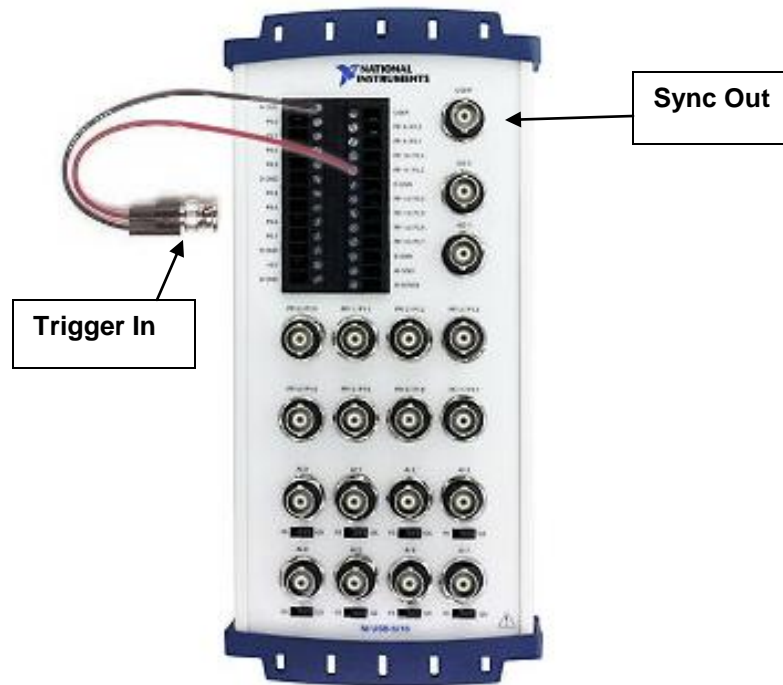


3. Wire a jumper from **PFI 12 / P2.4** to **USER** as shown below.

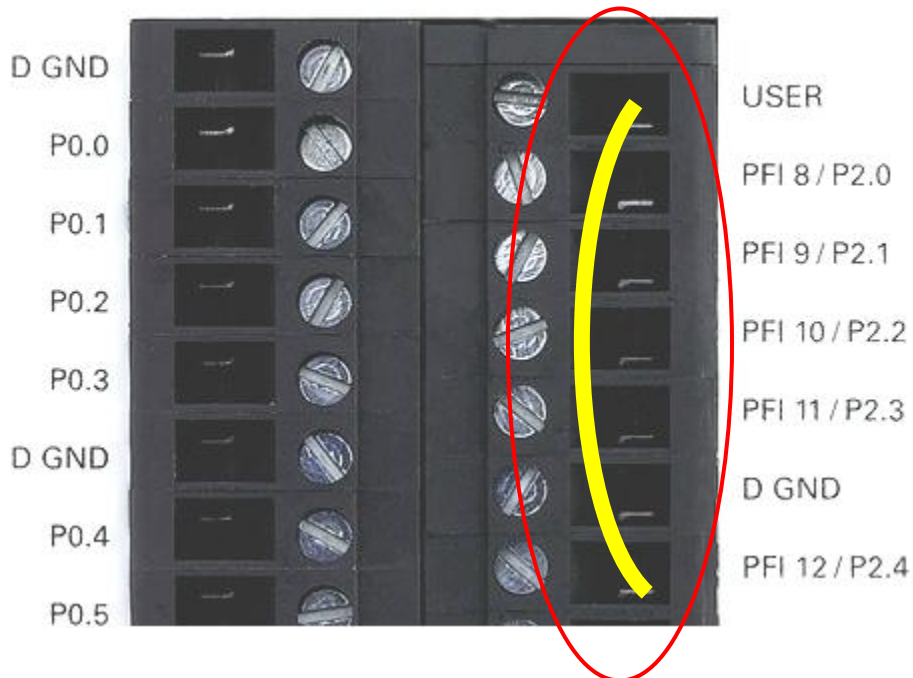

4. Wire a jumper from **P0.5** to **PFI 11 / P2.3**, as shown below.

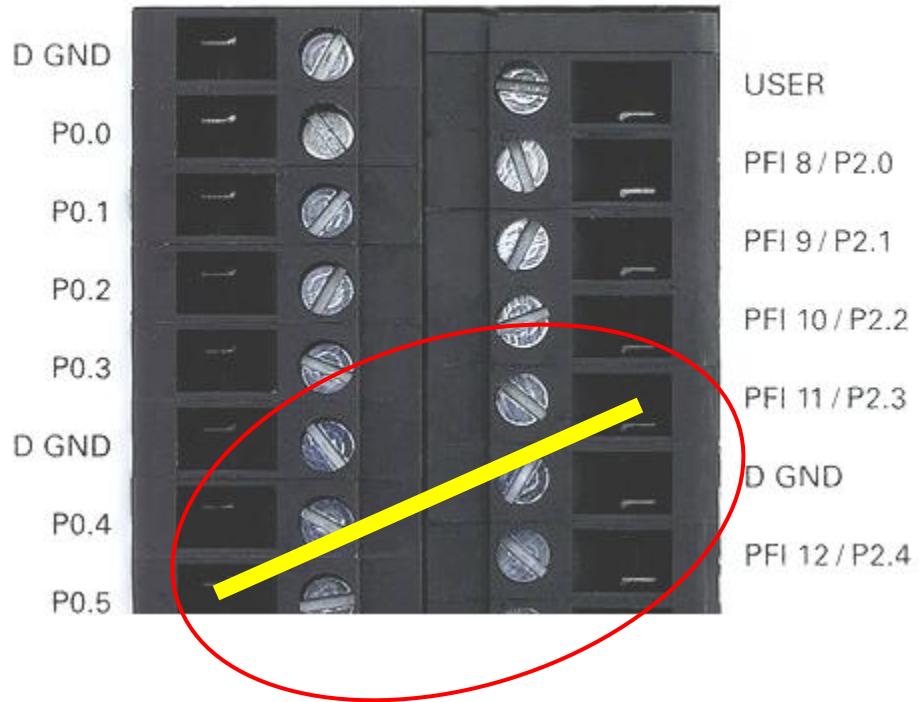

5. The picture below illustrates the wiring for the two jumpers.

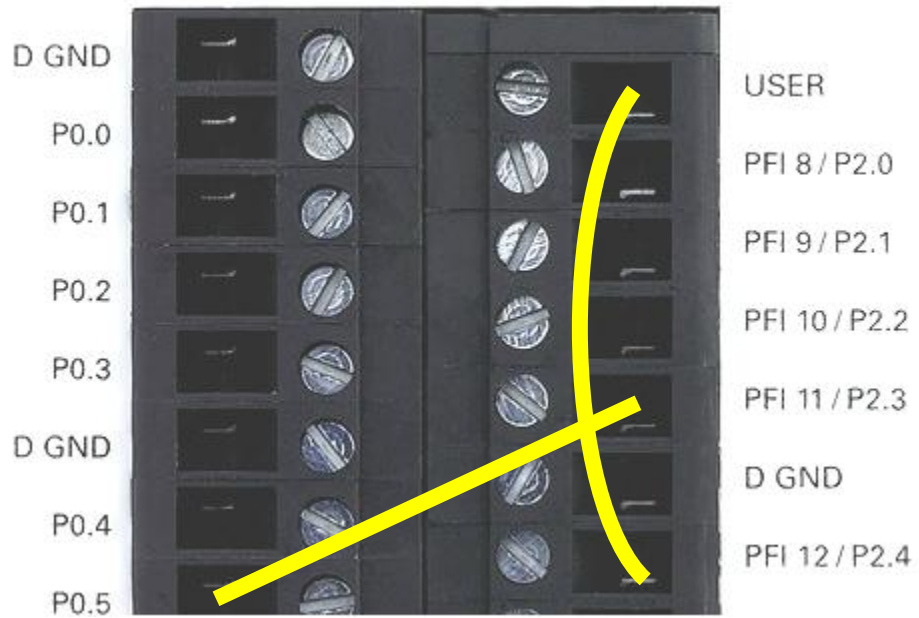

6. To start using your configured MiDAS DA hardware, please consult your *MiDAS DA User Guide*.

This application note is copyrighted by Xcitex Inc. and is supplied without specific warranty to any purpose and based on information currently available at the time of this writing. All specifications stated herein are subject to change without notice.

For further information on Xcitex products, visit www.xcitex.com or send an email to info@xcitex.com.

Xcitex Inc.
25 First Street, Suite 105
Cambridge, MA 02141 USA

