

# MIDAS DA

Version 1.4.0.1 – Release Notes

## General Description

---

This version update includes general stability and ease-of-use improvements. Additionally, we are proud to unveil a couple of exciting new features.

## Detailed Changes

---

1. Introduced the new Guided Demo Mode, which demonstrates recording and playback with step-by-step examples.
2. Introduced the new Data Summary Window.
3. Added Autoscaling option to Data Recording Windows.
4. Added support for Windows 7 x64.
5. Upgraded the National Instruments libraries.
6. Fixed bugs with Alignment Legend display bounds and initial state.
7. Fixed Samples/Frame changes not updating the Record Panel.
8. Fixed splash screen resolution conflicts.

## Installation Instructions

---

Run the installation executable and follow the on-screen instructions. The installation checks for Microsoft's .NET Framework v 2.0 as well as National Instrument's support for the .NET Framework and SafeNet's HASP software.

## Operating Systems

---

This version has been tested on Windows XP Professional Edition and Windows 7.

## **Hardware and Camera Compatibility**

---

MiDAS DA needs only an external trigger input and a sync. out interface to any camera and works easily with different brands. MiDAS DA supports E-Series and M-Series data cards from National Instruments and all accessories from National Instruments. Certain jumpers are required for proper operation.

## **Known Issues**

---

Portions of the electronic help file may not be up to date with respect to all recent changes.

## **How to get Technical Support**

---

Technical support information can be found on the Xcitex website on the Services and Support page at <http://www.xcitex.com/html/techsupport.php>. Contact information is available on the Contact Us page at <http://www.xcitex.com/html/contactus.php>. You may also contact Xcitex by phone, mail, or e-mail:

Xcitex, Incorporated  
25 First Street  
Suite 105  
Cambridge, MA 02141  
USA

US and Canada: 800-780-7836  
Telephone: 617-225-0080  
Facsimile: 617-225-2529  
[techsupport@xcitex.com](mailto:techsupport@xcitex.com)

Please have your key serial number, registration number and software version number available when calling.